
Water Sampling INSTRUCTIONS



Included Materials: Sampling bottles, sample ID stickers (two per water sample), chain of custody form, plastic bag (to keep form dry).

- Only test fixtures that are used for drinking, cooking, or making formula. Identify these water sources.
- Give each water fixture a unique name for easy identification. Write that name on the sample bottle AND the chain of custody form. Each sample has two numbered stickers. Place one sticker directly on the bottle and the other on the chain of custody form next to the fixture's name.

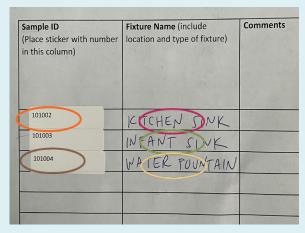



Collect COLD water samples from fixtures that have not been used for 8-18 hours. Collect first-draw samples in the morning after the water has been sitting still the night before. Collect the water sample immediately after turning on the faucet or valve. Do not allow any water to spill. Fill each sample bottle up to the 250ml mark on the bottle.



Sampling tips:

- **Do not** use the water fixture for 8-18 hours before taking the samples.
 - O To prevent people from accidentally using the water before sampling, you can tape off the area and post "do not use" signs.
- **Do not** use the facility's restrooms, sinks, or other water-using appliances or fixtures the morning before sampling as this can disrupt the water we want to sample.
- Water fixtures HAVE to be used regularly for accurate test results. Do not
 collect a sample from a fixture that has not been used in over 18 hours. If your
 facility is closed on the weekend, do not sample on Monday, and do not sample
 on the first day after a holiday or facility closure.
- **Do not** remove aerators or clean them before sampling.
- Complete the chain of custody form. Make a copy or take a photo for your records.

Water Shipping INSTRUCTIONS



Included Materials: Pre-filled FedEx shipping return label, box for shipping (use the same box your supplies arrived in), packing paper.

Place tightly sealed bottles in the box. Use the packing paper to keep the water samples from moving around.






Make sure the chain of custody form is complete and place in plastic bag.
Place the bag on top of sample bottles in box.

This form must be completed and returned for the lab to test your samples.



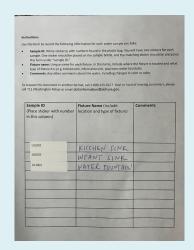

Remove old mailing labels. Seal box shut with packing tape, attach the pre-filled FedEx return label on box, and take box to the nearest FedEx location or call FedEx to have them come pick up the package: 800-463-3339 www.FedEx.com
Keep a copy of the tracking number. Shipping labels expire after 90 days.





- Send us an email at <u>leadfreekids@doh.wa.gov</u> so that we can notify the lab that samples are the way.
- We will e-mail your water test results. This typically takes 2-3 weeks after the lab receives your samples. Results will also be posted on our website quarterly.

Questions? Need Help? Contact <u>leadfreekids@doh.wa.gov</u> or 564-669-1098.